

How to download your NMR data to your computer.

Go to <http://www.versiontracker.com/dyn/moreinfo/macosx/15693> and download Fugu.

Open the dmg file, and copy Fugu to your applications folder.

Open Fugu.

On the menu bar, click Fugu -> preferences...

A window should open titled "Fugu Preferences: General"

Fill in the following information.

SFTP Server: chem9063.ucdavis.edu

Username: nmr

Initial Remote Directory: /home/archive/

Click "Set Defaults" and close the window.

Now click Connect, and when prompted, enter the password "1GoChem".

You should now be connected to the nmr server. There are 5 magnets listed in the archive, vnmrj300, vnmrj400, and vnmrj600 are the magnets in the basement, and nmr3470 is the magnet upstairs.

To transfer files, just locate the file or folder you want to download, and drag it from the window on the right to the window on the left.

Your newly created nmr data might not be available immediately, as the archive is backed up once an hour. For more information, check this sheet:

http://www.nmr.ucdavis.edu/NUTS_info/File_transfer_information.pdf or contact either John Berg or Jeff de Ropp.